AS.R ANALYTICAL BALANCES















Removable glass parts: side, top and back!

Quick access to information

Direct access to functions and databases is possible from the level of keyboard.

Database - a direct access to databasis Function - a direct access to the basic functions F1 to F4 - programmable function and navigation keys on the menu

ALIBI memory

The used ALIBI memory is a data secure area and allows to record up to 100 000 weighment records. It ensures security of constant data register in the long time period.

The AS.R series represents a new standard level for analytical

Additionally, the new R series balances by means of pictograms signal the activated working mode, connection with the Internet, the battery charge level, balance service functions. Also a number of displayed measuring units has been increased.

The balance precision and the measurement accuracy is assured by automatic internal adjustment, which takes into consideration temperature changes and time flow.

AS.R series balances feature several communication interfaces: 2 x RS 232, type A USB, type B USB and optional WiFi. The housing is made of plastic, and the pan is made of stainless steel.

DATABASES IN R SERIES BALANCES

In new AS.R series balances the information system is based on 5 databases, which allows for several users to work with several products databases, and the registered weighing results can be subject to further analysis.

The data is registered in 5 databases:

- users (up to 10 users),
- products (up to 1000 products),
- weighments (up to 5000 weighments),
- -tares (up to 100 tares),
- -ALIBI memory (up to 100 000 weighments).

There is two directions data exchange within the system thanks to a quick USB interface. New balances allow to import and export databases using USB pen drives.

R series balances fulfill GLP requirements.

balances. They feature a new, readable LCD display which allows a clearer presentation of the weighing result. Besides, the display has a new text information line allowing to show additional messages and data, e.g. product name or tare value.



Parts counting



Dosing



Checkweighing



Percent deviations



Statistics



Animal weighing



Pipettes calibration

Control



Statistical Quality



Autotest (GLP, Filter)



GLP procedures



Under hook weiahina



Totalizing



Density determination



Peak hold



Packaged Goods Control



Newton unit measurement



Replaceable



Technical specification: AS 62.R2 AS 60/220.R2 AS 82/220.R2 M M Max capacity 62 g 60 q / 220 q 82 q / 220 q Minimum load 1 mg 1 mg 1 mg Readability 0,01 mg 0,01 mg / 0,1 mg 0,01 mg / 0,1 mg -220 g Tare range -62 g -220 g $0,015 \text{ mg } (Rt \le 2 \text{ g})$ $0,015 \text{ mg } (Rt \le 2 \text{ g})$ $0,015 \text{ mg } (Rt \le 2 \text{ g})$ $0.02 \text{ mg} (2 \text{ g} < \text{Rt} \le 50 \text{ g})$ $0.02 \text{ mg} (2 \text{ g} < \text{Rt} \le 50 \text{ g})$ Repeatability * $0.02 \text{ mg} (2 \text{ g} < \text{Rt} \le 50 \text{ g})$ $0.03 \text{ mg} (50 \text{ g} < \text{Rt} \le 60 \text{ g})$ $0.03 \text{ mg} (50 \text{ g} < \text{Rt} \le 82 \text{ g})$ $0.03 \text{ mg} (50 \text{ g} < \text{Rt} \le 60 \text{ g})$ $0,1 \text{ mg } (60 \text{ g} < \text{Rt} \le 220 \text{ g})$ $0.1 \text{ mg } (82 \text{ g} < \text{Rt} \le 220 \text{ g})$ \pm 0,06 mg (to 60 g) ± 0,06 mg (to 82 g) Linearity ± 0,06 mg $\pm 0.2 \text{ mg} (60 \text{ g} \div 220 \text{ g})$ ±0,2 mg (82 g ÷ 220 g) Pan size open-work pan Ø 90 mm or (Ø 85 mm - option) Working temperature +10° ÷ +40°C Relative air humidity ** 40% ÷ 80% Stabilization time 6 s / 3.5 s 6 s $6 \, s / 3,5 \, s$ Sensitivity drift 1 ppm/°C in temperature +10° ÷ +40°C Minimum weight (USP) 30 mg 30 mg 30 mg Minimum weight (U = 1%, k = 2) 3 mg 3 mg 3 mg Interface 2 × RS 232, USB-A, USB-B, WiFi - option Power supply*** 12 ÷ 16 V DC / 2,1 A Adjustment/calibration internal (automatic) Display LCD (backlit) Net weight/Gross weight 5,4 kg / 7,5 kg 5,4 kg / 7,5 kg 5,4 kg / 7,5 kg Packaging size 495 × 400 × 515 mm

^{*} Repeatability is expressed as a standard deviation from 10 weighing cycles.


^{**} Non-condensing conditions


^{*** 250} mA for balances without WiFi module, 350 mA for balances with installed WiFi module

Technical specifica	ation:			
	AS 110.R2	AS 160.R2	AS 220.R2	AS 310.R2
	M	M	M	M
Max capacity	110 g	160 g	220 g	310 g
Minimum load	10 mg	10 mg	10 mg	10 mg
Readability	0,1 mg	0,1 mg	0,1 mg	0,1 mg
Tare range	-110 g	-160 g	-220 g	-310 g
Repeatability *	0,1 mg (Rt ≤ 110g)	0,1 mg (Rt ≤ 160g)	0,1 mg (Rt ≤ 220g)	0,1 mg (Rt ≤ 220g) 0,2 mg (220g < Rt ≤ 310g
Linearity	± 0,2 mg	± 0,2 mg	± 0,2 mg	± 0,3 mg
Pan size	Ø 100 mm	Ø 100 mm	Ø 100 mm	Ø 100 mm
Working temperature	+10° ÷ +40°C			
Relative air humidity **	40% ÷ 80%			
Stabilization time	3,5 s			
Sensitivity drift	1 ppm/°C in temperature +10° ÷ +40°C			
Interface	2 × RS 232, USB-A, USB-B, WiFi - option			
Power supply***	12 ÷ 16 V DC / 2,1 A			
Adjustment/calibration	internal (automatic)			
Display	LCD (backlit)			
Net weight/Gross weight	5,4 kg / 7,5 kg	5,4 kg / 7,5 kg	5,4 kg / 7,5 kg	5,4 kg / 7,5 kg
Packaging size	495×400×515 mm			

^{*} Repeatability is expressed as a standard deviation from 10 weighing cycles.

Dimensions:

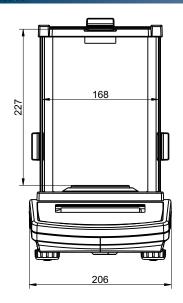

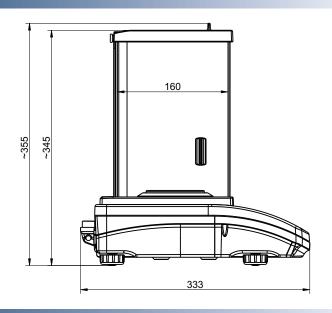

Accessories:

Antivibration table SAL/STONE	Density determination kit	
Rack for under hook weighing	Additional LCD display "WD-6"	
Professional weighing table	Power adapter with battery and charger ZR-02	
Impact printer Epson	PC keyboard USB	
Printer USB PCL	External USB memory (FAT files format)	
Holders for glass vessels	Mass standard	
"PW-WIN" computer software	Adjustment weight	
"RAD-KEY" computer software	USB A- USB B cable (balance - computer, balance - PLC printer)	
Antistatic ionizer DJ-02	Cable RS 232 (scale - computer) "P0108"	
Bar code scanner	Cable RS 232 (scale, Epson, Citizen printer) "P0151"	
Bar code scanner USB HID	"Tare" or "Print" foot button	



^{**} Non-condensing conditions

 $^{^{\}star\star\star}$ 250 mA for balances without WiFi module, 350 mA for balances with installed WiFi module